

Flow Solution™ FS 3700 Automated Chemistry Analyzer

Available Cyanide by Ligand Exchange
OIA-1677-09 by Flow Injection Analysis (FIA)
Cartridge Part Number 330092CT

Scope and Application

This method is used for determining available cyanide in water and wastewater by ligand exchange, flow injection analysis, and amperometric detection according to **USEPA Method OIA-1677-09** (Reference 1, 5, 6, 9) and **ASTM Method D 6888-09** (Reference 2). This method is used in the USEPA's data gathering and monitoring programs associated with the Clean Water Act, Resource Conservation and Recovery Act, Comprehensive Environmental Response, Compensation and Liability Act, and Safe Drinking Water Act.

This method is also applicable to **ASTM D 7284-08** Total Cyanide by manual distillation gas diffusion amperometry and **USEPA OIA 1677-DW** "free" cyanide.


Cyanide ion (CN⁻), hydrogen cyanide in water (HCN(aq)), and cyano-complexes of zinc, copper, cadmium, mercury, nickel, and silver can be determined by this method. The presence of polysulfides and colloidal material may prove intractable for application of this method.

Method Performance

Range	2.0 µg/L–5.0 mg/L
Rate	30 samples/hour
Precision	3% RSD
Method Detection Limit (MDL)	0.5 µg/L

The range may be extended to analyze other concentrations by changing the size of the sample loop. The quality of the analysis is assured through reproducible calibration and testing of the Flow Injection Analysis (FIA) system.

A general flow diagram of the Flow Injection Analysis (FIA) system is shown below (see Figure 1 for a detailed flow diagram).

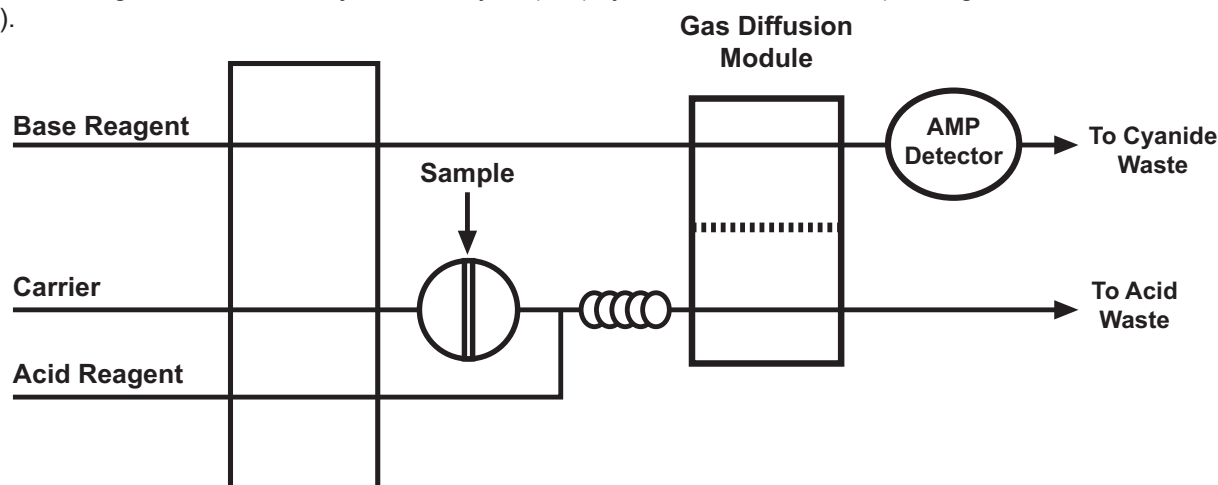

Reagents and Calibrants

Chemical Name	CAS #	Chemical Formula	Part Number
Acetic acid, glacial	64-19-7	$C_2H_4O_2$	
Acetone	67-64-1	C_3H_6O	
Bismuth nitrate pentahydrate	10035-06-0	$Bi(NO_3)_3 \cdot 5H_2O$	
Water, deionized		H_2O	
5-[4-(Dimethylamino) benzylidene] rhodanine	536-17-4	$C_{12}H_{12}N_2OS_2$	
DOWFAX 2A1	12626-49-2		A000080
Ethylenediamine	107-15-3	$C_2H_8N_2$	
Ligand exchange reagent, available cyanide (WAD) reagent A			A001416
Ligand exchange reagent, available cyanide (WAD) reagent B			A001417
Mercury (II) cyanide	594-04-1	$Hg(CN)_2$	
Nickel potassium cyanide	14220-17-8	$K_2Ni(CN)_4$	
Potassium cyanide	151-50-8	KCN	
Silver nitrate	7761-58-3	$AgNO_3$	
Sodium acetate	127-09-3	$C_2H_3O_2Na$	
Sodium hydroxide	1310-73-2	NaOH	
Sulfide carrier reagent (SCR)			A001668
Sulfuric acid, concentrated	7664-93-9	H_2SO_4	
Sulfuric acidification reagent (SAR)			A001872
WAD carrier			A001125
WAD AR			A001501

Summary of OIA-1677-09

OIA-1677-09 summarizes the method and interferences as follows below (Reference 6).

Method

- Prior to analysis, treat the sample to remove potential interferences. Add ligand exchange reagents to the sample. Thermodynamically stable complexes form with the transition metal ions listed in, releasing the cyanide ion from the cyano-complexes. Inject an aliquot of the treated sample into the FIA system (Reference 2). Addition of acid converts the cyanide ion to hydrogen cyanide gas (HCN), which passes under a gas diffusion membrane. The hydrogen cyanide gas diffuses through the membrane into an alkaline receiving solution where it converts back to cyanide ion. The cyanide ion is monitored amperometrically with a silver working electrode, silver/silver chloride reference electrode, and platinum/stainless steel counter electrode at an applied potential of zero volt. The current generated is proportional to the cyanide concentration present in the original sample.